For staining polyacrylamide gels

## Silver Stain KANTO III



What kind of reagents have you been using for staining polyacrylamide gels? Please try "Silver Stain Kanto III" which is a simple and fast kit for silver staining.

## Fast staining

It takes only one hour to stain gels after electrophoresis, which is faster compared with conventional methods.

## High reproducibility

This kit provides the clearer staining image with high reproducibility.

# Re-use after CBB staining

Gels stained with CBB can be re-stained using this kit.

## High sensitivity

The sensitivity for detection is 10-100 fold for proteins and 5-10 fold for nucleic acids compared with CBB method and Ethidium Bromide method, respectively.

### Ready-to-use

This kit is ready-to-use only by dilution of each solution with de-ionized water.

## Safety

Since sodium hydroxide or ammonia water is not used, there is no danger to generate explosive silver amide.

### **♦**Components**♦**

Reagent	Principal component	content
A (Stock solution for amplification)	Dithiothreitol	200ml × 1
B (Stock solution I for staining)	Silver nitrate	200ml × 1
C (Stock solution II for staining)	Silicotungstic acid	200ml × 1
D (Stock solution I for development)	Formaldehyde	200ml × 1
E (Stock solution II for development)	Potassium sodium carbonate	200ml × 1
F (Termination solution)	Citric acid	200ml × 1

#### **♦**Procedure**♦**

#### **Solution for fixation:**

Mix methanol 50ml, acetic acid 10ml and de-ionized water 40ml.

#### Solution of amplification:

Mix Reagent A 5ml and de-ionized water 95ml.

#### **Solution for staining:**

Mix Reagent B 5ml, Reagent C 5ml and de-ionized water 90ml.

#### Solution for development:

Mix Reagent D 5ml, Reagent E 5ml and de-ionized water 90ml.

#### **Solution for termination:**

Reagent F is to be used as it is.

STEP	TIME
fixation	20 min
amplification	10 min
washing	10 min
staining	10 min
washing	1 min
washing	1 min
development	5-10 min
termination	10 min

#### **♦**Product**♦**

Cat. No.	Name	Size
37937-96	Silver Stain KANTO III	1 Kit (For 80x80mm 40gels)

### igspaceRelated productsigspace

Cat. No.	Name	Size
36535-96	Rapid CBB KANTO	1L





試薬部

(03) 3663-7631 541-0048 大阪市中央区瓦町2丁目5番1号 (06) 6231-1672 812-0007 福岡市博多区東比恵2丁目22番3号 (092) 414-9361

<< http://www.kanto.co.jp e-mail: reag-info@gms.kanto.co.jp >>