

CicaAllerTest[®] House dust Mite allergen Der f 1



You can detect
house dust mite allergen
Easily!

For Der f 1




Dermatophagoides farinae and *Dermatophagoides pteronyssinus* are recognized as the most important indoor environmental allergens worldwide. From these mites, various allergen proteins have been identified. Among them, the allergen (Der 1) contained in stool is recognized as a universal standard of the degree of mite pollution and has a significant correlation with the number of mites. CicaAllerTest[®] House dust Mite allergen Der f 1 is an immunochromatography reagent that qualitatively detects Der f 1, allergen from stool of *D. farinae*.

For Identification of contaminated areas of *D. farinae*.
Confirmation of cleaning condition.
Evaluation of the performance of antiallergenic agents.


Features

- **Short time!** - Decision time is only 15 minutes.
- **High sensitivity!** - Minimum detection sensitivity is 1 ng / mL (Der f 1).
- **Simple!** - No special equipment is required.


Operation method

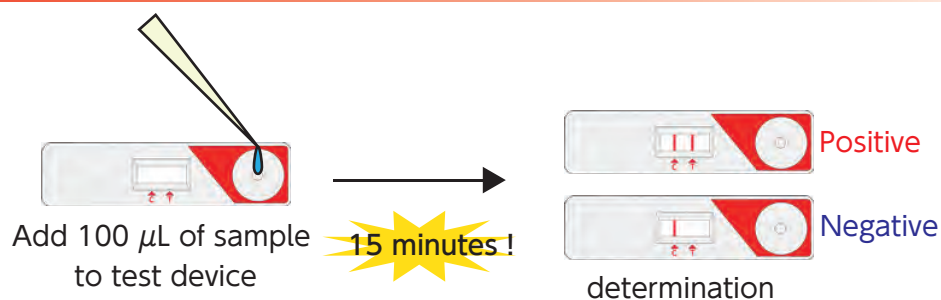

Determination method

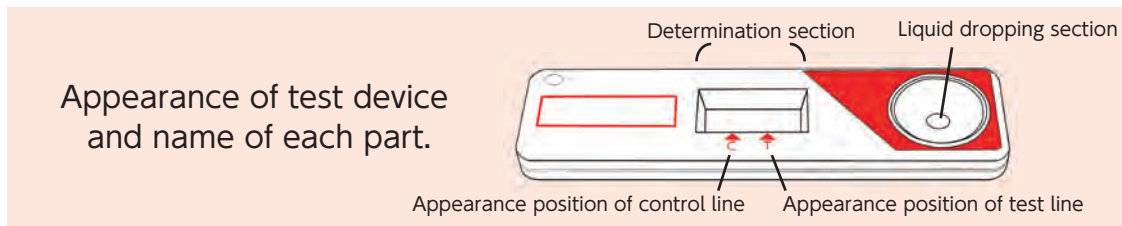


Result is judged by the presence or absence of the control line and the test line of the determination section.

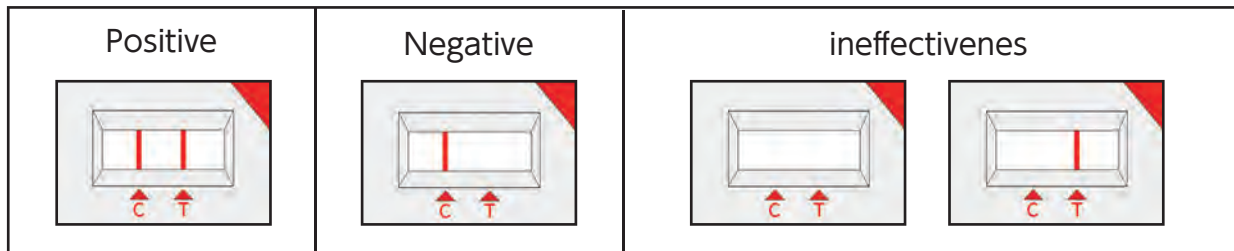

- ※ Please keep decision time (15 minutes) strictly.
- ※ This reagent is for qualitative measurement, it is not developed for quantitative purposes. In order to quantify, please use other methods such as ELISA.

Examples of test

Different concentration of Der f 1 was tested.

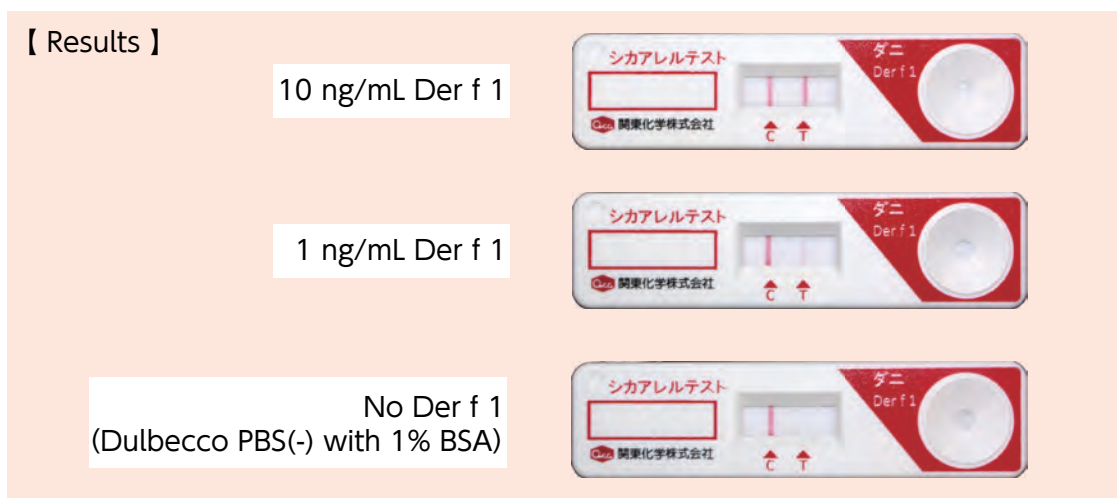

Product Information

Product No.	Product Name	Package size
07961-67	CicaAllerTest® House dust Mite allergen Der f 1	10 tests

KANTO CHEMICAL CO., INC.
REAGENT DIVISION

East Muromachi Mitsui BLDG, 2-1, Nihonbashi Muromachi 2-chome,
Chuo-ku, Tokyo, 103-0022, JAPAN

Telephone +81-3-6214-1092

Telefax +81-3-3241-1053

<http://www.kanto.co.jp>

E-mail: kanto-61@gms.kanto.co.jp