

BMP7 selective inhibitor

| | |
|----------------------|-------------|
| Code | PG-005 |
| MW (Acetate): | 2085.29 |
| Purity: | ≥95% (HPLC) |
| Amount: | 5 µg |

Formulation: This vial contains 5 µg of the titled compound, which has been lyophilized from 0.2 µm filtered DMSO solution.


Reconstitution: Reconstitute at 30~100 µM in DMSO. Centrifuge the vial before opening.

<Example>

50 µM solution prepared when 5 µg peptide reconstituted by 48 µL DMSO.

Storage: Store at -20°C or below for the container vial unopened. Recommend to use the peptide as quickly as possible after opening and reconstitution.

Activity: Measured by its ability to inhibit smad1 phosphorylation induced by BMP4 or BMP7 in A549 cell.

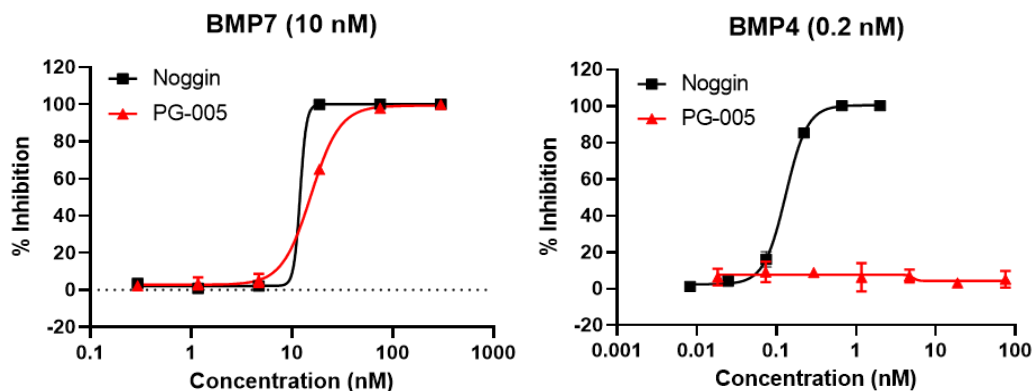

This product is "For Research Use Only"